
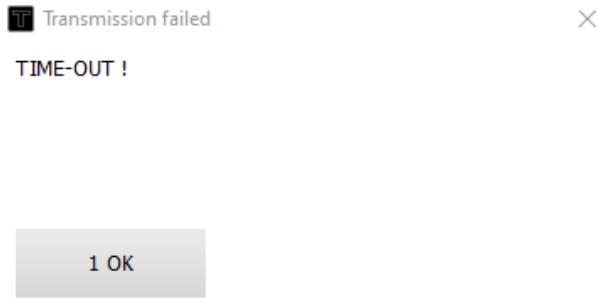
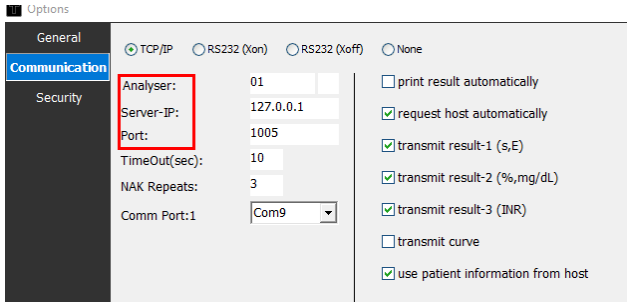


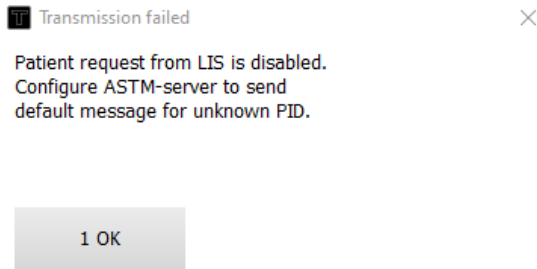

Malfunction TECAM does not connect to ASTM Server

Problem-1: 	Solution: Configure TECAM to correct IP,port of LIS-Server 
Problem-2: 	Solution: TECAM sends a test request ('Q') for PID=123456789000 to LIS-Server. The LIS must reply with a default profile Add a default order for PID= "123456789000" into your LIS or configure the LIS to send a default profile , if PID is unknown. 
===== TECAM send test request 'Q' for PID=123456789000 <pre>TECO <ENQ> LIS <ACK> TECO <STX>1H \^& 01^ P 97 20250122120053<CR><ETX>ED<CR><Lf> LIS <ACK> TECO: <STX>2Q 1 ^123456789000<CR><ETX>87<CR><Lf> LIS <ACK> TECA <STX>3L 1 N<CR><ETX>06<CR><Lf> LIS <ACK> TECO <EOT></pre>	===== LIS sends test order 'O' for PID=123456789000 <pre>LIS [ENQ] TECO [ACK] LIS [STX]1H \^& 01^ [CR][ETX]18 [CR][LF] TECO [ACK] LIS [STX]2P 1 1 Inf1^^^ 19641231 [CR][ETX]2A [CR][LF] TECO [ACK] LIS [STX]3O 1 123456789000 ^1^1^3 R [CR][ETX]E2 [CR][LF] TECO [ACK] LIS [STX]4L 1 N [CR][ETX]07 [CR][LF] TECO [ACK] LIS [EOT]</pre>